

LabTemp® 20/40

STORAGE / INOCULATION / INCUBATION



STAGE 1 STORAGE

Upon receipt, place LabTemp® 20 on agar plate (agar side up). Store at 2-8°C (in the dark).

LabTemp® 20/40 packaged ready for use




STAGE 2 INOCULATION

Allow the medium to warm to room temperature then inoculate (LabTemp® 20 will indicate when to inoculate 17-20°C)

STAGE 3 INCUBATION

Place inoculated agar medium with LabTemp® 40 (LabTemp® 40 on the agar side up) into incubator set at 35°C ± 2.0°C in aerobic atmosphere for 18-48 hours. When removing plate from incubator check LabTemp® 40 to see if agar plate is at proper temperature.

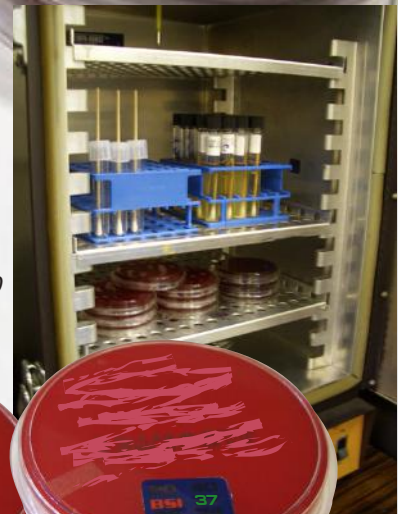

After sufficient (temperature/time) incubation, plate shows isolated colonies of organisms.



BIOSYNERGY, INC.

TIME / TEMPERATURE / SOLUTIONS

1940 E. DEVON • Elk Grove Village, IL 60007
800-255-5274 OR 847-956-0471